


Pre-grown, ready-to-use biofilm test carriers
for antimicrobial performance testing

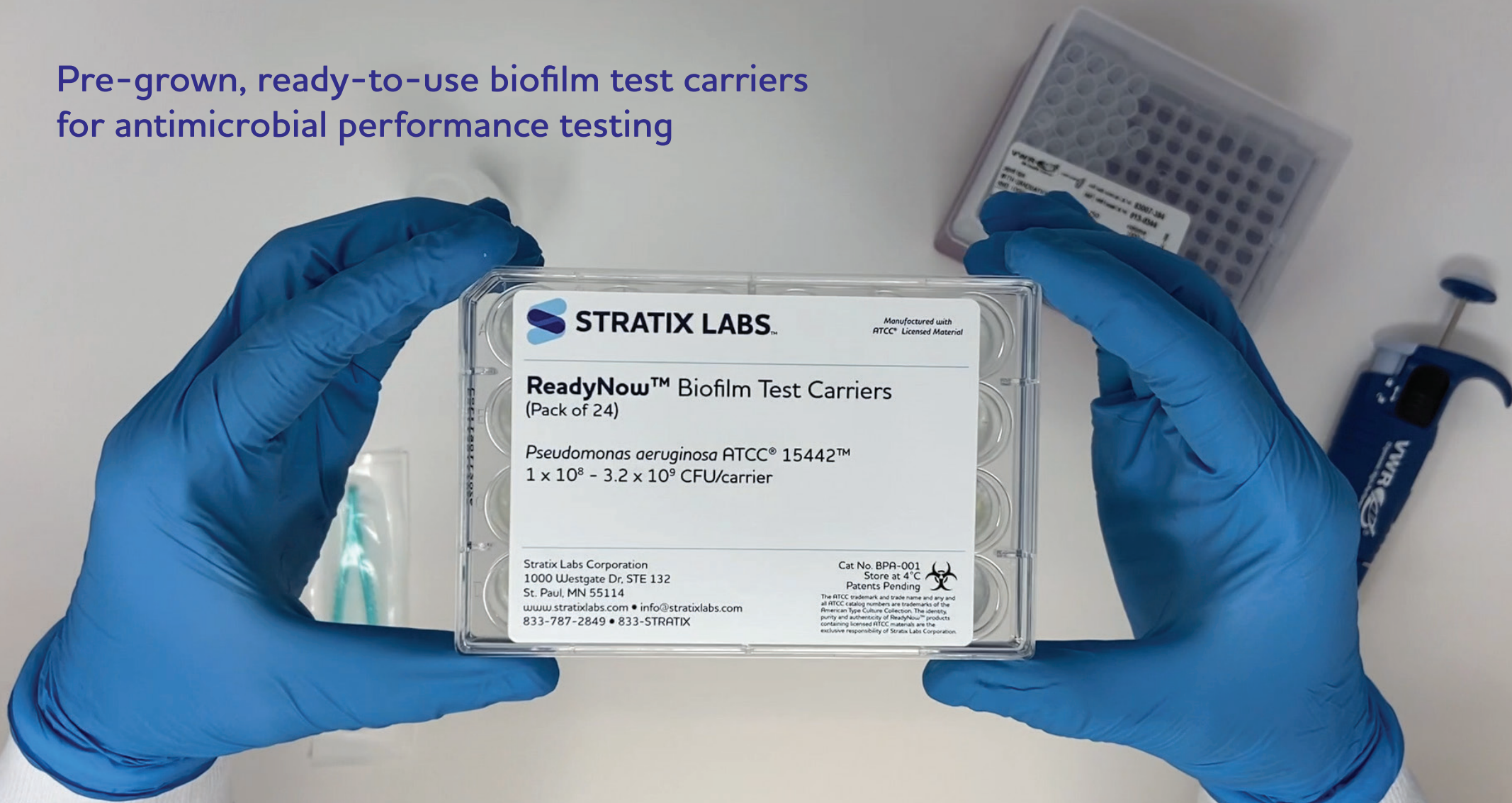

Manufactured by:

Stratix Labs Corporation • 1000 Westgate Dr, STE 132 • St. Paul, MN 55114
www.stratixlabs.com • sales@stratixlabs.com • Tel: (833) 787 2849; (833) STRATIX

Distributed in partnership with:

BioSurface Technologies Corporation • 421 W Griffin Drive • Bozeman, MT 59715
www.biofilms.biz • sales@biofilms.biz • Tel: (866) 306-2620



*We take care of the prep work,
so you don't have to.*



ReadyNow™ Biofilm Test Carriers

No reactors. No pumps. Ready for testing. Simply rehydrate for 15 minutes at room temperature, and you're ready to test for anti-biofilm efficacy.

Each kit includes 24 pre-grown, ready-to-use biofilm test carriers and a 50 mL bottle of rehydration solution.

Available Organisms

Pseudomonas aeruginosa ATCC® 15442™
1.0 x 10⁸ - 3.2 x 10⁹ CFU per carrier

Staphylococcus aureus ATCC® 6538™
1.0 x 10⁷ - 3.2 x 10⁸ CFU per carrier

Manufactured by:

Stratix Labs Corporation • 1000 Westgate Dr, STE 132 • St. Paul, MN 55114
www.stratixlabs.com • sales@stratixlabs.com • Tel: (833) 787 2849; (833) STRATIX

Distributed in partnership with:

BioSurface Technologies Corporation • 421 W Griffin Drive • Bozeman, MT 59715
www.biofilms.biz • sales@biofilms.biz • Tel: (866) 306-2620




***Ready-to-use in less than 20 minutes.
Anti-biofilm efficacy testing has never been easier.***

1




Rehydrate

2

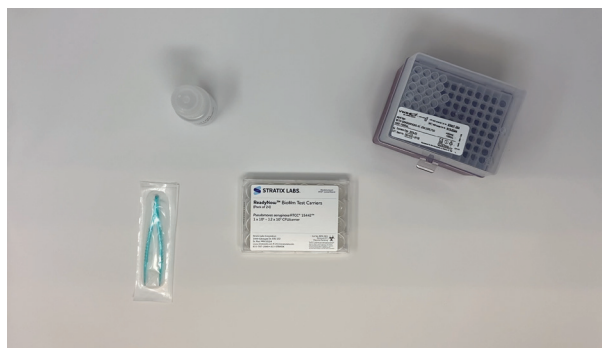

Wait 15 minutes

3

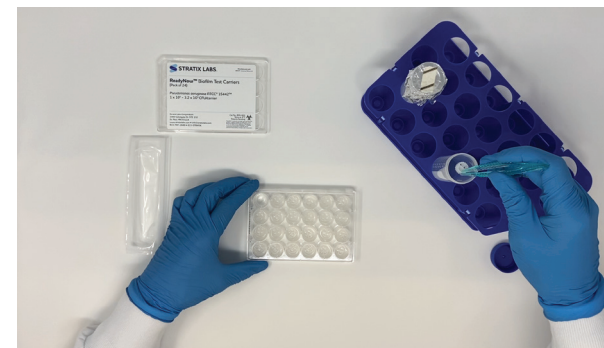


Conduct your assay



Precise & In-spec Mean Log CFU/carrier: 8.54


Pseudomonas aeruginosa ATCC® 15442™

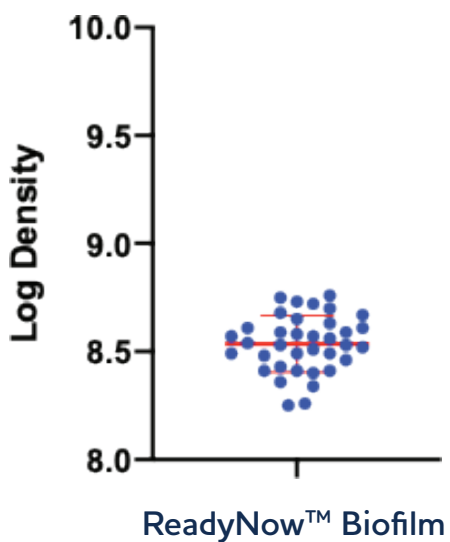

Mean	8.54
Std dev.	0.13
Max.	8.76
Min.	8.25
n	36

Responsive & Repeatable Disinfectant Efficacy Test

Pseudomonas aeruginosa ATCC® 15442™

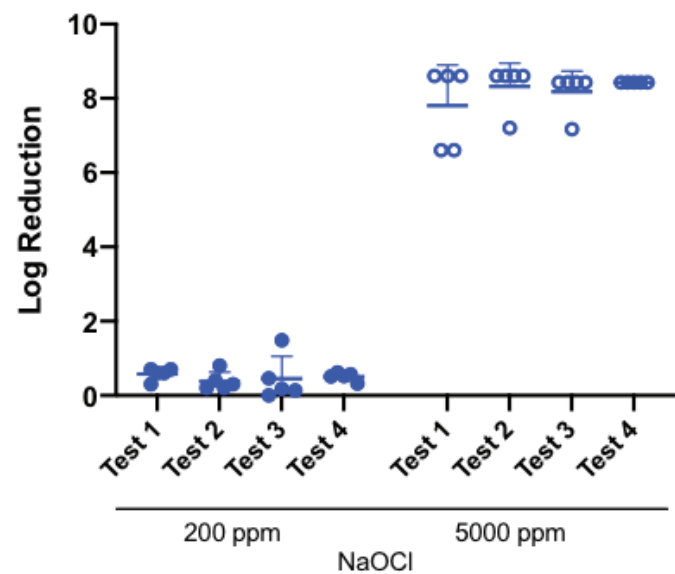


Disinfectant Efficacy Testing						
NaOCl						
Test Day	200 ppm			5000 ppm		
	Mean	SD	n	Mean	SD	n
1	0.58	0.16	5	7.80	1.10	5
2	0.38	0.24	5	8.32	0.63	5
3	0.45	0.60	5	8.18	0.56	5
4	0.51	0.11	5	8.43	0.00	5
Total	0.48	0.32	20	8.18	0.68	20



Comparison of Time Required to Prepare Biofilm Test Samples ReadyNow™ Biofilm Test Carriers vs Manual Preparation with Biofilm Reactor



98% LESS
PREP TIME

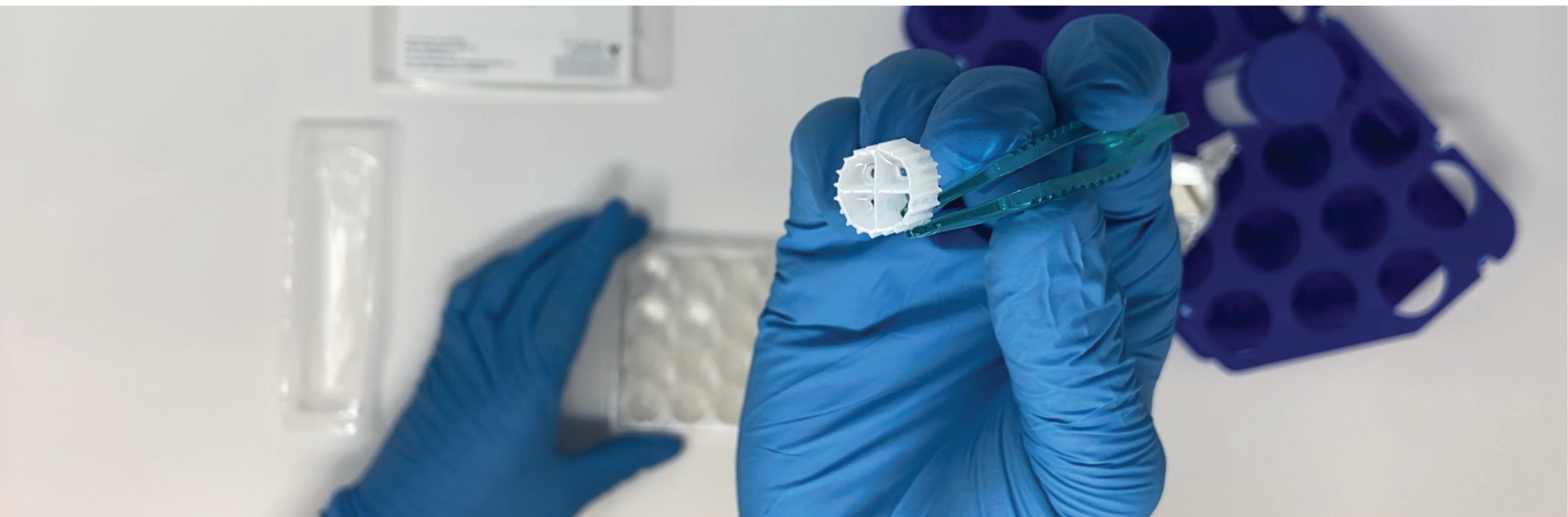

Manufactured by:

Stratix Labs Corporation • 1000 Westgate Dr, STE 132 • St. Paul, MN 55114
www.stratixlabs.com • sales@stratixlabs.com • Tel: (833) 787 2849; (833) STRATIX

Distributed in partnership with:

BioSurface Technologies Corporation • 421 W Griffin Drive • Bozeman, MT 59715
www.biofilms.biz • sales@biofilms.biz • Tel: (866) 306-2620